

# pulseCheck NX

## The Modular Autocorrelator

Pulse Measurement Perfection  
with the Multitalent from APE

The pulseCheck platform grows with your tasks. Suitable for characterizing virtually all ultrafast pulse lasers, APE's pulseCheck autocorrelator covers a wide range of wavelengths and pulse widths. This flexibility is achieved by using interchangeable optics sets, typically consisting of a nonlinear crystal and a dedicated detector module.



- Exchangeable Optics Sets for wavelength coverage from 200 nm to 12  $\mu\text{m}$
- Wide range of sensitivity levels covered with PMT, PD, and TPA
- High temporal resolution with a sampling resolution down to 50 attoseconds
- Toggle between interferometric and intensity autocorrelation
- Integrated Controller for data acquisition, data fitting, and fast FPGA data processing
- Software aided wavelength tuning of crystal (phase matching)
- Fully compliant with DIN 58175-2:2021-04 (pulse measurement)
- NX-Software and USB as well as Ethernet connection
- TCP/IP remote control with standardized command set for easy programming
- Option: FROG for complete pulse characterization


# pulseCheck NX Overview

## Overview

Specifications	pulseCheck NX 50	pulseCheck NX 150	pulseCheck SM 2000
Pulse Width	5 fs - 15 ps	10 fs - 40 ps	20fs - 500 ps
Wavelengths	200 nm - 12 $\mu\text{m}$ , depending on optics set	200 nm - 12 $\mu\text{m}$ , depending on optics set	200 nm - 12 $\mu\text{m}$ , depending on optics set
Delay Resolution	50 attoseconds	200 attoseconds	1 femtosecond
Refresh Rate	10 Hz	7.5 Hz	Depending on scan range, 120 ps/sec
Intensity Resolution	18 bit	18 bit	18 bit
Detector Types	PMT, TPA, PD	PMT, TPA, PD	PMT, TPA, PD
Electronics	Integrated	Integrated	Integrated
Software	NX Software	NX Software	V 2019 Software
Connectivity	USB, Ethernet	USB, Ethernet	USB, Ethernet
			see page 14 pulsecheck SM 2000 for details

## High Resolution & Strong Processing Performance

The advanced system-on-chip based architecture enables measurements with low latency, a high time resolution down to 50 attoseconds and an intensity dynamic range of 18 bit. This allows full access to even the tiniest autocorrelation features.

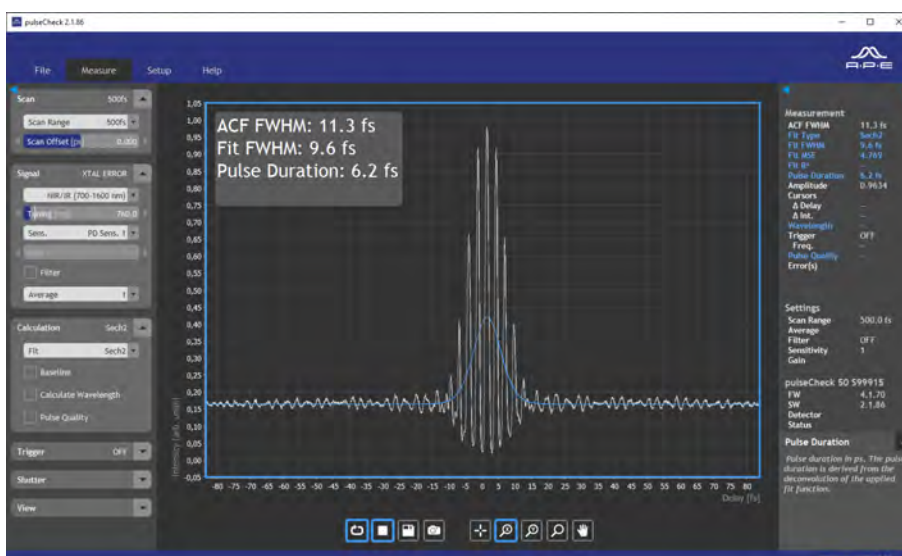

Software GUI for pulseCheck NX


## ... Flexibility for your Experiments

### Maximum Functionality through Modular Design

APE meets the growing need for maximum functionality and flexibility with the modular concept on which the pulseCheck autocorrelator series is based. Thus, Optics sets to upgrade the wavelength range or FROG capability can be added any time.

### From Ultrashort to Longer Pulses

Various pulseCheck configurations can be adapted to the individual needs of pulse duration measurement. The measurement of extra long pulse durations up to 500 ps is possible with pulseCheck SM.

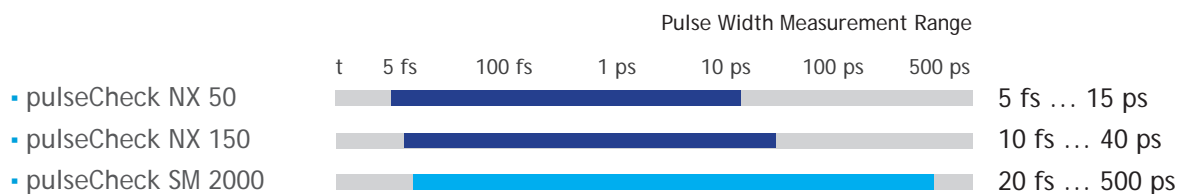

### High Sensitivity and Low Noise with Three Types of Detectors

- Photodiode Detector (PD)
- Photomultiplier (PMT)
- Two-Photon Absorption (TPA)

Standard sensitivity up to  $1 \text{ W}^2$   
 Highest sensitivity up to  $10^{-6} \text{ W}^2$   
 High sensitivity up to  $10^{-2} \text{ W}^2$




# pulseCheck NX

## Optics Sets

A variety of different and exchangeable optics sets cover a wide wavelength range: from UV at 200 nm to mid IR at 12  $\mu\text{m}$ .


## Photodiode (PD)

- **Typ. Sensitivity:**  $1 \text{ W}^2$
- **Rep. Rate:**  $> 10 \text{ Hz}$
- **Measurement Mode:** Collinear and Non-collinear

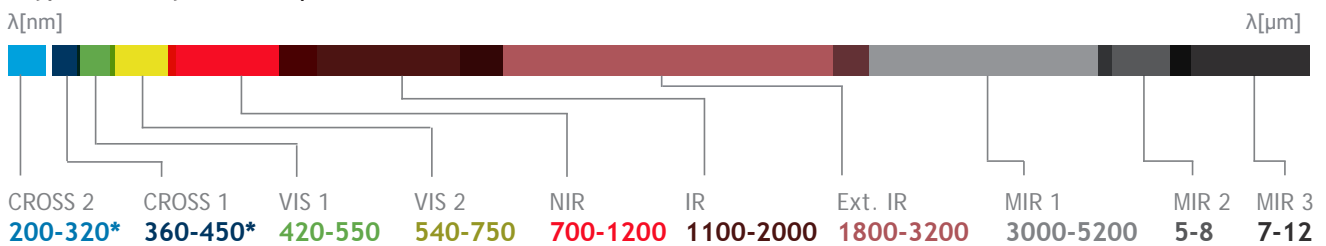


## Photomultiplier (PMT)

- **Typ. Sensitivity:** up to  $10^{-6} \text{ W}^2$
- **Rep. Rate:**  $> 250 \text{ kHz}$
- **Measurement Mode:** Collinear and Non-collinear

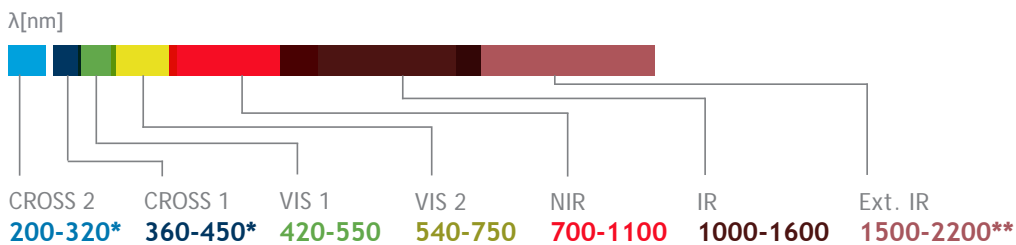

## Two Photon Absorption (TPA)

- **Typ. Sensitivity:**  $< 0.1 \text{ W}^2$  (NIR/IR) |  $< 1 \text{ W}^2$  (VIS) |  $< 100 \text{ W}^2$  (UV)
- **Rep. Rate:**  $> 10 \text{ Hz}$
- **Measurement Mode:** Collinear Intensity

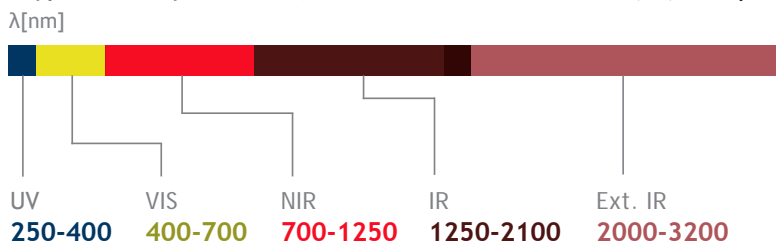

\* For cross-correlation: wavelength range depends on pump wavelength

\*\* For the wavelength range 1500 - 2200 nm we recommend to use the highly sensitive IR detector "Extended IR PD SELECTED"

# pulseCheck NX Specifications

## Specifications

Measurable Pulse Width Range*	5 fs - 15 ps with pulseCheck NX 50; 10 fs - 40 ps with pulseCheck NX 150
Wavelength Range	200 nm - 12 $\mu$ m, depending on Optics Set
Exchangeable Detector (Optics Sets)	PMT, PD, or TPA
Dealy Sampling Resolution	Down to 50 attoseconds (NX 50), 200 attoseconds (NX 150)
Refresh Rate	10 Hz (NX 50); 7.5 Hz (NX 150)
Sensitivity	Typically $1 \dots 10^{-6} W^2$ depending on Optics Set**
Intensity Resolution	18 Bit
Pulse Repetition Rate	PD, TPA: 10 Hz and above; PMT: 250 kHz and above
Type of Measurement Mode	PMT, PD : non-collinear intensity, collinear interferometric; TPA: hybrid non-collinear intensity, collinear interferometric;
SHG Tuning for Phase Matching	PMT/PD: automatic; TPA: not applicable
Trigger Mode	TTL for $f < 50$ kHz
Input Polarization	Linear horizontal, vertical available as option
Input Beam Coupling	Free-space (6 mm aperture); Option: fiber coupling FC/PC, FC/APC, SMA
Max Input Power	1 W (e.g. oscillator with a rep. rate of approx. 70 MHz) or 10 $\mu$ J (e.g. amplified system with rep. rates in the kHz range), whichever results in lower value
Connectivity	USB and Ethernet
Remote Control	Possible via TCP/IP (SCPI command set)
Calibration	NIST traceable calibration certificate included
Calibration	NIST traceable calibration certificate included

## NX Software & Electronics

Electronics	<ul style="list-style-type: none"> <li>• Integrated controller for data acquisition, processing and data fitting</li> <li>• Fast FPGA data post-processing</li> <li>• &gt; 10k datapoints per scan</li> <li>• Connectivity: USB and Ethernet</li> <li>• In-field software and firmware updates</li> </ul>
Software	<ul style="list-style-type: none"> <li>• Intuitive and easy-to-navigate</li> <li>• Plug &amp; play installation</li> <li>• Real-time display of the pulse width (measured ACF and fitted pulse width)</li> <li>• Fitting Routines: Gaussian, Sech<sup>2</sup>, Lorentz</li> <li>• Measurement of central wavelength</li> </ul>
NX Software Enhancements	<ul style="list-style-type: none"> <li>• The metadata dashboard displays all measurement results and device status</li> <li>• The new help section displays hints and explanation for all software buttons and functions</li> <li>• The new monochromatic Contrast Screen Mode helps to enhance the readability of the screen while wearing laser safety glasses</li> </ul>
Remote	<ul style="list-style-type: none"> <li>• Direct remote control via TCP/IP (no PC software required)</li> <li>• Users can create their own measurement remote routines using the standardized SCPI command set and data acquisition</li> </ul>

\* 13 fs<sup>2</sup> internal dispersion at 800 nm for short-pulse option to be considered

\*\* Sensitivity is defined as measured sensitivity including Optics Set, defined as average power times peak power of the incident pulses  
 $P_{AV} * P_{peak}$

## Options

---

- Various Optics Sets
- Fiber coupling
- Polarization rotator
- FROG

## Compliance

---

- CE certified
- NIST traceable calibration with HeNe laser
- Fully compliant with DIN 58175-2:2021-04 (Autocorrelation measurement methods)

## Dimensions and Power

---

Dimensions	244 x 201 x 309 mm (pulseCheck NX 50) 244 x 201 x 395 mm (pulseCheck NX 150) See appendix for technical drawings
------------	--

Power	95 ... 240 V, 50 ... 60 Hz, 60 W
-------	----------------------------------

## pulseCheck SM 2000 Long-Range Autocorrelator

Pulse Width Measurement  
up to 500 ps

With the revised version of the long-range delay technology in pulseCheck SM, the new autocorrelator offers a scanning range up to three times larger and a measuring speed up to five times faster than the previous long-range models. The pulseCheck SM includes a high performance controller that is integrated into the device.

The pulseCheck SM is suitable for particularly long pulse durations of up to 500 ps. At the other end, short pulses down to 20 fs can be measured. A resolution of 1 fs and a sampling rate of 1 MHz is available over the entire measurement range.



- Pulse widths from 20 fs up to 500 ps
- Optics Sets from 200 nm to 12  $\mu\text{m}$
- Measurement speed of 120 ps/sec
- Automatic phase matching
- Integrated high performance controller
- Toggle between interferometric and intensity autocorrelation
- Wide range of sensitivity levels covered with PMT, PD, and TPA
- USB & Ethernet connectivity and TCP/IP remote control
- Gaussian,  $\text{Sech}^2$ , and Lorentzian fitting routines
- NIST traceable calibration

# pulseCheck SM 2000 Specifications

## Specifications

Measurable Pulse Width*	20 fs ... 500 ps (20 fs - 150 ps directly measured, 150 ps - 500ps half ACF measured)
Wavelength Range	200 nm - 12 µm, depends on Optics Set
Optics Sets	Exchangeable; Detector-Types: PMT, PD, or TPA
Delay Resolution	1 fs
Delay Linearity	< 1 %
Scan Speed	120 ps/s
Sampling Rate	1 MHz
Sensitivity**	Typically $1 \dots 10^{-6} \text{ W}^2$ depending on Optics Set
Recommended Repetition Rate	PD, TPA: 1 Hz ... GHz; PMT: 250 kHz ... GHz
Type of Measurement Mode	PMT, PD : non-collinear intensity, collinear interferometric; TPA: hybrid collinear intensity
Trigger Mode	TTL, repetition rate 1 Hz - 10 kHz
Input Polarization	Linear horizontal, vertical available as option
Input Beam Coupling	Free-space (6 mm aperture); Option: fiber coupling FC/PC, FC/APC, SMA
Max Input Power	1 W (e.g. oscillator with a rep. rate of approx. 70 MHz) or 10 µJ (e.g. amplified system with rep. rates in the kHz range), whichever results in lower value
Software	Included; Real-time display of data and pulse parameters
Fitting Routines	Gaussian, $\text{Sech}^2$ , Lorentz
Connection	USB, Ethernet
Remote Control	Possible via TCP/IP (SCPI command set)
Calibration	NIST traceable calibration certificate included

## Dimensions and Power

Dimensions	253 x 190 x 572 mm (W/H/L) See appendix for technical drawings
Power	95 ... 240 V, 50 ... 60 Hz, Power 60 W

\* 13 fs<sup>2</sup> internal dispersion at 800 nm for short-pulse option to be considered

\*\* Measured sensitivity including Optics Set, defined as average power times peak power of the incident pulses  $P_{AV} * P_{peak}$




# Appendix Technical Drawings

All Dimensions in mm

## pulseCheck NX

- Multitalent for any task

page 3

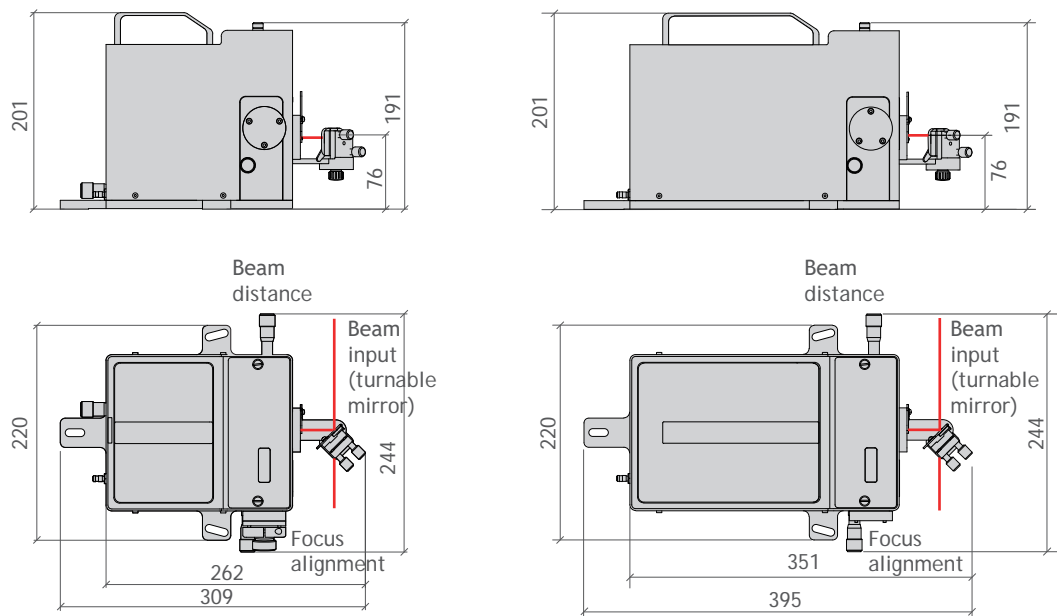


## pulseCheck SM 2000

- Long-range autocorrelator

page 9

**PHOTO TECHNICA** [www.phototechnica.co.jp](http://www.phototechnica.co.jp)  
**フォトテクニカ株式会社**  
 〒336-0017 埼玉県さいたま市南区南浦和 1-2-17  
 TEL:048-871-0067 FAX:048-871-0068  
 e-mail:voc@phototechnica.co.jp

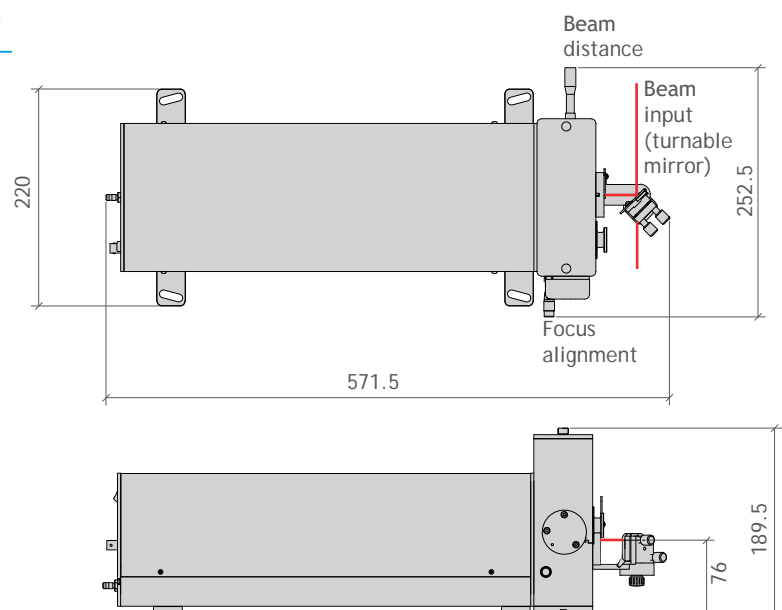

### Contact

APE Angewandte Physik & Elektronik GmbH  
 Plauener Str. 163-165 | Haus N | 13053 Berlin | Germany  
 T: +49 30 986 011-30 F: +49 30 986 011-333  
 E: sales@ape-berlin.de www.ape-berlin.de